



LAUDA Service-Info ECO cooling thermostats

Warning 044 – ‘Chiller missing’

	<p><u>Safety information:</u></p> <p>Repair and service work should be done by qualified personnel only:</p> <ul style="list-style-type: none">• Service on electronics must be done by a trained electrician.• All refrigeration and cooling service must be done by a trained and certified refrigeration technician. <p>Disposal of the used refrigerant should be done by a trained and certified refrigeration technician only!</p> <p>Remove mains plug before opening the device!</p>
	<p>This manual should only be regarded as a collection of helpful solutions to technical problems concerning LAUDA thermostats and cooling units.</p> <p>No responsibility is taken for the correctness of this information.</p>

Symptom: After short operation in cooling mode the warning message 044 ‘Chiller missing’ is displayed.

Problem: Missing cable connection between device underpart and control head or bad connection due to not fully inserted plug into socket 51H.

Solution: Check if the plug of the connection is inserted properly.
Please follow the next step if it looks like on picture 1:

Switch off device and push the plug with some more force whilst counter with the other hand at the front panel. Switch on the device again.

